

VDA 278 Analysis using Thermal Desorption with a Dynatherm-GC/MS System

The German Automobile Association method VDA 278 is a thermal desorption analysis for organic emissions from interior components of a vehicle. This analysis deals with a VOC value (up to C20-Eicosane) based on Toluene, and a FOG value (C16-C32) based on C16 (Hexadecane). It deals with the emissions coming from non-metallic molded parts occurring in motor vehicles. Materials such as textiles, adhesives, foam, leather, and similar substances are included. This method is used for both qualitative, as well as, semiquantitative evaluation of the degassed emissions. The VOC emissions are reported semiquantitatively as toluene, whereas the FOG value is reported as condensable substances found in the materials as hexadecane. Single substances of the emission can also be identified and quantitated using this method or a variation of this method.

A serial dilution was made of a Kovats VOC standard (100µg/ml) to give a final concentration of 10ng/µl in CS₂. A thermal desorption tube (6") packed with Tenax was spiked with the standard VOC mix (20ng/2µl). The tube was thermally desorbed for 280°C/30min on the Dynatherm 9300 TDA interfaced to a GC/MS (Figure 1 and Figure 3). GC parameters were established to resolve C8 (octane) to C20 (Eicosane) and C8 to C 27 (for FOG analysis). A corresponding analysis of carpet fibers was performed by placing them into an empty 6" desorption tube and desorbing them for 30 minutes at 90°C (VOC portion, Figure 2) and for 60 minutes at 120°C (FOG portion, Figure 4). A number of aliphatic hydrocarbons were identified (C13,C14,and C15) in the VOC por-

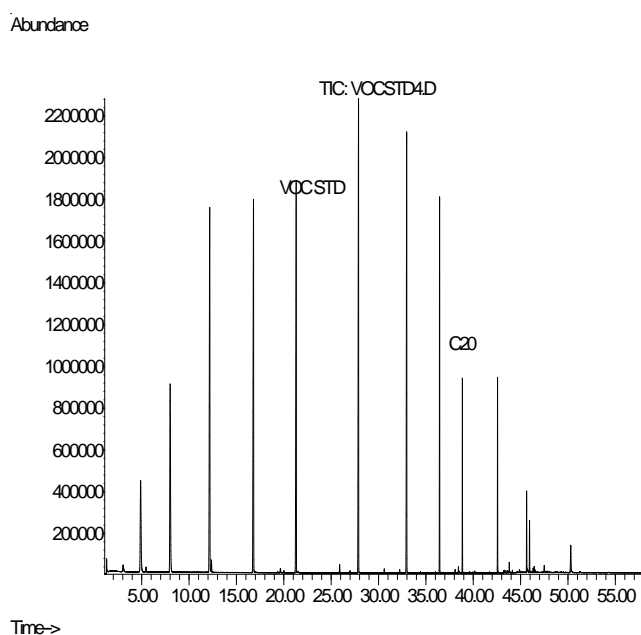


Figure 1

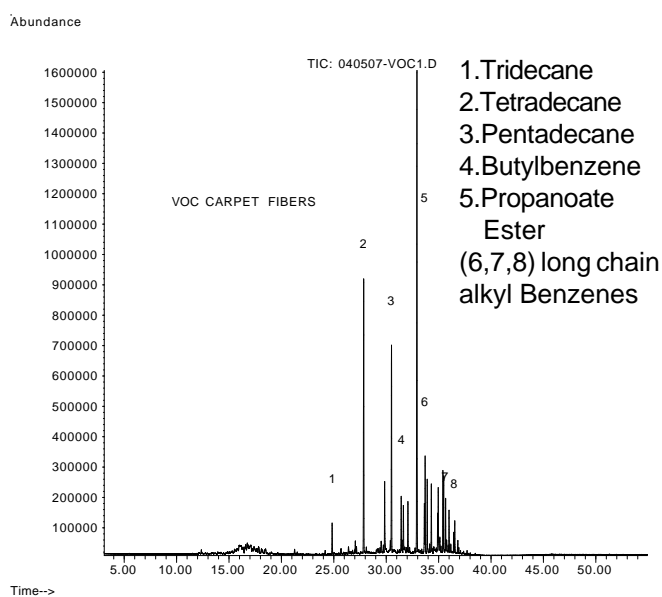


Figure 2

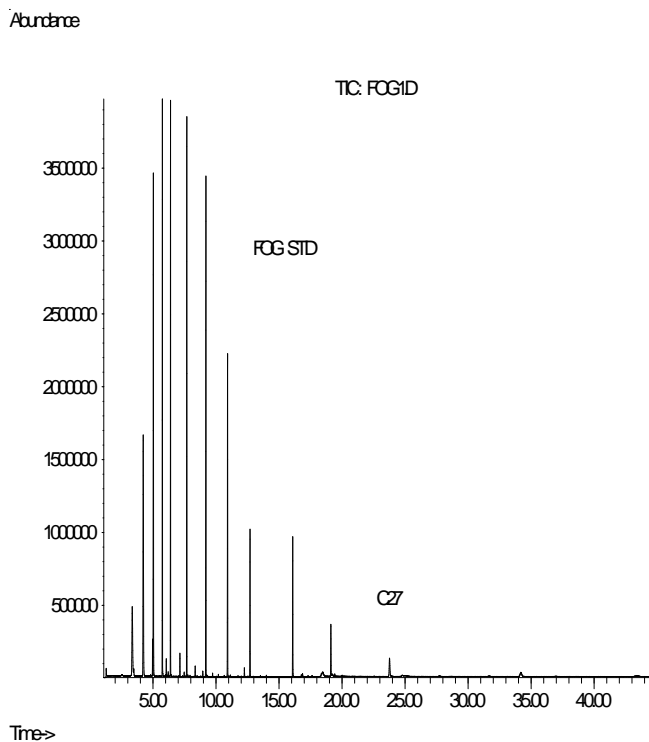


Figure 3

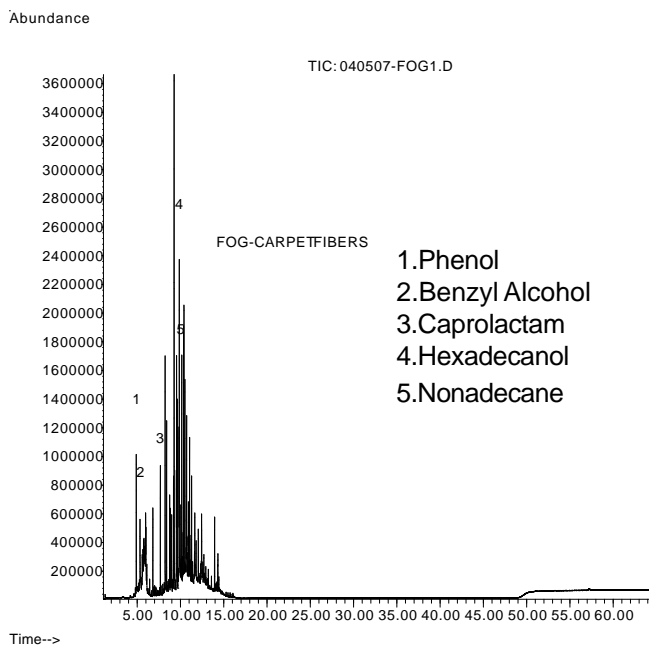


Figure 4

1. Phenol
2. Benzyl Alcohol
3. Caprolactam
4. Hexadecanol
5. Nonadecane

tion as well as a propanoate ester and a number of long chain alkyl benzenes. The FOG fraction had a number of aromatics, phenol, benzyl alcohol, as well as, caprolactam. There were also long chain aliphatic alcohols and several alkanes detected.

GC/MS Method:

VOC Portion:

Initial: 40°C/2min
 Rate: 3°C/min to 92°C, 5°C/min to 160°C, 10°C/min to 280°C, hold for 10min

FOG Portion:

Initial: 50°C/2min
 Rate: 25°C/min to 160°C, 10°C/min to 280°C, hold for 30min

TDA 9300 Conditions:

Kovats Standard:

Valve Oven: 280°C
 Transfer Line: 280°C
 Tube Heat: 280°C/30min
 Trap Heat: 280°C/10min

VOC (Carpet Fibers) TDA 9300 Conditions:

Valve Oven: 280°C
 Transfer Line: 280°C
 Tube Heat: 90°C/30min
 Trap Heat: 280°C/20min

FOG (Carpet Fibers) TDA 9300 Conditions:

Valve Oven: 280°C
 Transfer Line: 280°C
 Tube Heat: 120°C/60min
 Trap Heat: 280°C/20min

CDS Analytical, Inc. has been a leader in the design and manufacture of laboratory instruments for sample preparation and analysis since 1969. We are dedicated to providing the best possible instruments for both research and routine analysis. Well known in the field of pyrolysis, CDS manufactures the Pyroprobe® 5000, 5150, 5200 and 5250 autosampler for the introduction and analysis of solid materials by GC, MS and FT-IR. CDS offers a complete line of dynamic headspace instruments for the analysis of volatile organic compounds in environmental, pharmaceutical and food applications, including the model 8400 four-position autosampler for complex, multicomponent materials investigation. Our customers, their requirements and applications are important to us. To help meet your needs, we offer a wide range of analytical information and the services of our applications laboratory. If you would like additional information, please contact us at the address below, call us at 1 800 541 6593, or log onto www.cdsanalytical.com.